


ML230 Focused-ultrasonicator


Uninstallation and Packaging Instructions for Return Shipment

Instrument Shutdown Procedures

1. After decontamination of the instrument (see decontamination protocol), ensure the following:
 - a. Water bath is completely drained and has been blotted dry (using soft, lint-free cloth or KIM wipes).
 - b. Waste bottle is empty and has been thoroughly rinsed.
 - c. Supply bottle is empty and has been thoroughly rinsed.
2. Close SonoLab by touching the **"White X"** at the upper right corner of the screen.
3. When the **"Drain Bath?"** pane appears, touch **"Exit SonoLab"**.
 - a. Water bath will have drained after following the decontamination procedure.

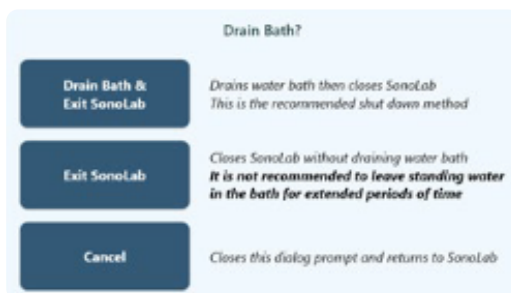

4. Once the **"Power Off Computer?"** pane appears, touch **"Exit SonoLab & Power Off Computer."**

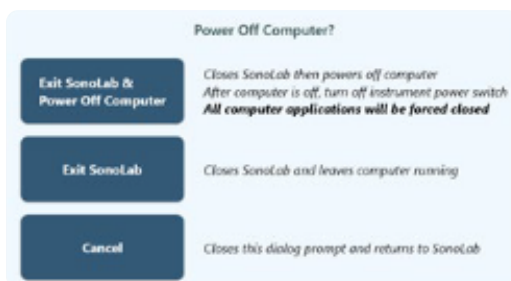


5. Once the computer is off, move the ML230 power switch to the **"OFF"** position.
 - a. Power switch is located on the rear panel of the ML230, about halfway up the left side, above the power cord.


Instrument Uninstallation Procedures

NOTE: When lifting the instrument, use both sides of the base to lift, DO NOT lift by the plastic parts at the front of the instrument.


1. Unplug the AC power cord from the power outlet and then from the ML230 power inlet (located about halfway up the left side of the ML230 rear panel).

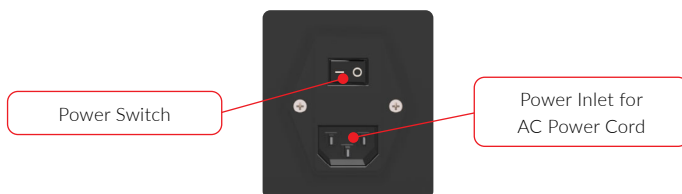


2. Remove the ML230 safety cover.
3. Unscrew the sample tray assembly's thumbscrew from the motion system and then slide out the sample tray.

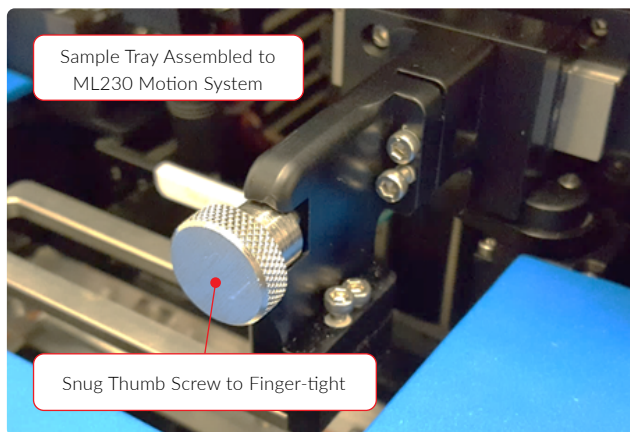

4. If the supply and waste bottles are still attached to the ML230, press the quick disconnect port and slide out of the water control bay.

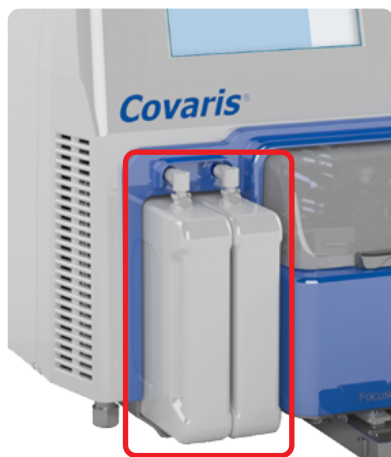

5. Before placing the instrument in the cardboard shipping box, tighten each leveling foot so it is snug and completely shortened.
- a. Shorten both legs on the right side of the instrument by turning clockwise.
 - b. Shorten both legs on the left side of the instrument by turning clockwise.

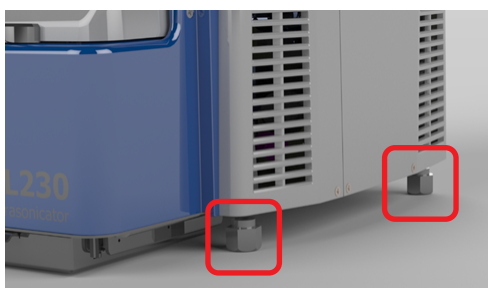

Instrument Packaging Procedures

1. Place the instrument in the foam base.



4. Place box around the instrument packaging assembly.




2. Place bag on the top of the instrument.



5. Place supply bottle, waste bottle, and all other accessories into the remaining accessory cardboard box and place above the instrument foam top.

6. Place the sample tray into its respective foam and place [preferably] inside the accessory box.

3. Place foam top on the instrument.

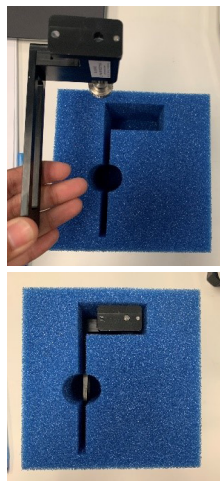

7. Place the lid to close off the assembly.
8. Place shipping label on the side of the box.

Notes

Contact Covaris

- Telephone
 - United States: Tel: +1 781.932.3959 during the hours of 9:00 a.m. to 5:00 p.m., Monday through Friday, Eastern Standard Time (EST), Greenwich Mean Time (GMT-05:00)
 - Europe: Tel: +44 (0) 845 872 0100, during the hours of 9:00 a.m. to 5:00 p.m., Monday through Friday, Greenwich Mean Time
- E-mail queries to applicationsupport@covaris.com or techsupport@covaris.com
- For questions regarding Demos, please contact customerservice@covaris.com (US) or eucustomerservice@covaris.com (EU).

Information subject to change without notice. For research only. Not for use in diagnostic procedures.

USA: Tel: +1 781.932.3959 | Email: customerservice@covaris.com | Europe: Tel: +44 (0)845 872 0100 | Email: emeacustomerservice@covaris.com
APAC: +86 137 6427 6714 | Email: APACcustomerservice.com | Web: www.covaris.com | Applications: applicationsupport@covaris.com
Service and Support: techsupport@covaris.com | 2022© Covaris, LLC