

96 Shuttle RFID Label Mounting Instructions

Trademarks

Patented or patent pending and registered or registration-pending trademarks of Covaris are protected.

Registered names and trademarks used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

WCS and AFA are either registered trademarks or trademarks of Covaris, LLC in the United States and/or other countries.

Information subject to change without notice.

For Research Use Only (RUO).

Not for use in Diagnostic Procedures.

Product covered by Patent No. US6, 719,449, US6, 948,843, and other pending applications.

2022© Covaris, LLC. All Rights Reserved.

Covaris, LLC

14 Gill St.

Woburn, Massachusetts

01801-1721 USA

Technical Support

Ongoing assistance with the operation or application of the equipment and/or troubleshooting is provided via:

1. Telephone:

- US: +1 781.932.3959 during the hours of 9:00 a.m. to 4:00 p.m., Monday through Friday, Eastern Standard Time (EST), Greenwich Mean Time (GMT-05:00)
- EU: +44 (0)845 872 0100, during the hours of 9:00 a.m. to 5:00 p.m., Monday through Friday, Greenwich Mean Time (GMT-05:00)


2. E-mail instrumentation queries to techsupport@covaris.com and application queries to applicationsupport@covaris.com

Table of Contents

96 Shuttle RFID Label Mounting Instructions	4
Revision History	5

96 Shuttle RFID Label Mounting Instructions

1. Remove the old RFID label by peeling it off the shuttle by grabbing the exposed edge of the RFID label and peeling sideways.
2. Install a new RFID label on the shuttle in the exact position shown below, in the exact orientation shown below.
3. Run finger along top face of RFID label over areas with metal behind it while applying light pressure to adhere label to shuttle.

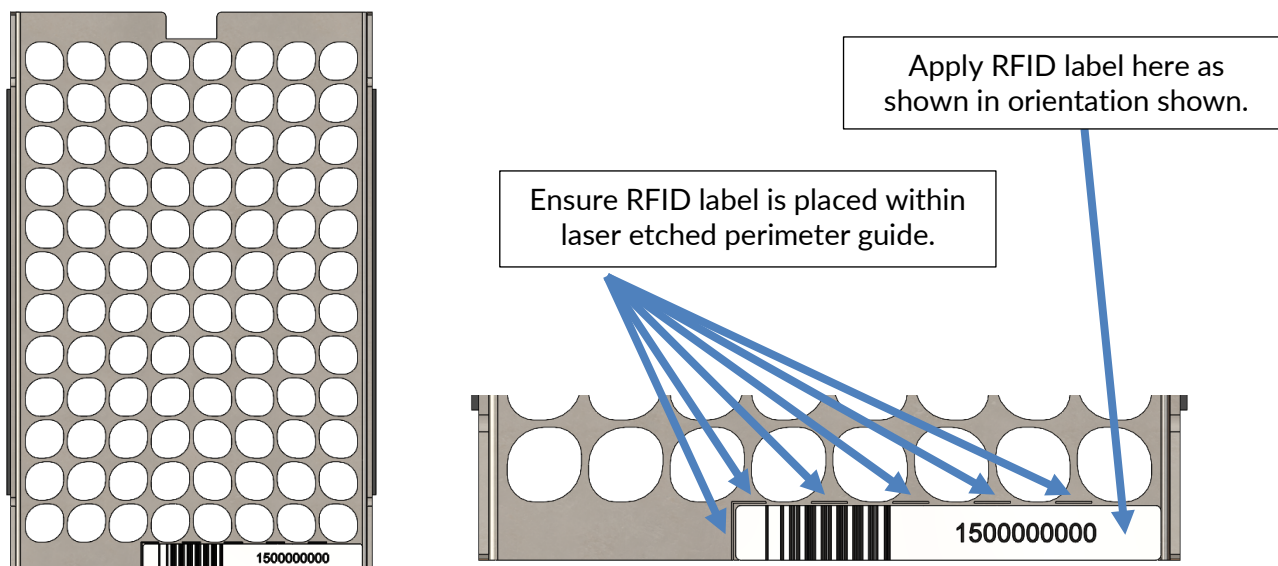

Installation example shown below. Always use the label with the RFID marker on it!



**Do not place label off of perimeter of shuttle or curl it upwards in any way.
It must lay flat along entire length of label. Precision is important.**

Notes

N/A

Revision History

Document Number	Revision	Date	Description of Change
010612	A	01/2023	Initial release of document